



Application Note

## Vitroids™ and Lenticule® discs

How to use certified reference materials (CRMs) to perform a spiking protocol for method verification.

## Case

Spiking protocol preparation to reach a low level of Salmonella inoculation:

3-5 CFU/test portion as requested in ISO 16140-3 (Protocol for the verification of reference methods and validated alternative methods) when applying Protocol #3.

Choose an appropriate CRMs, for this case VT000292 Salmonella enterica subsp. enterica serovar Abony WDCM 00029 is selected, as an example. This product has a mean value range between 15-80 CFU.

To verify which is the precise value of a specific lot, its <u>certificate</u> must be checked.

From the certificate of lot BCCG2046, as an example

Sample: Salmonella enterica subsp. enterica serovar Abony WDCM 00029 VT000292		
Certified value (geometric mean value)	Expanded uncertainty (log <sub>10</sub> value)	Expected range
4.9E+01 cfu per disc	0.03	2.4E+01 - 9.8E+01 cfu per disc

In this case the certified value is 49 CFU of Salmonella per disc.

## **Protocol**

- 1. Prepare the initial suspension of the sample to be inoculated (25 g sample + 225 mL diluent).
- 2. In a separate vial, rehydrate one vitroid in 10 mL Buffered Peptone Water (BPW). Do not spend more than 10-15 minutes for rehydration and spiking (point 3), because the bacteria can start growing in BPW. If the reference material contains 49 cfu per disc (with uncertainty: from 24 cfu to 98 cfu), this Salmonella suspension will contain 4,9 cfu/mL (with uncertainty from 2,4 to 9,8 cfu/mL).
- 3. Spike the initial suspension (25 g sample + 225 diluent) with 500 microliters of the homogenized Salmonella suspension. Considering the reference value, the sample finally would be spiked with a reference value of 2,45 cfu/25 g sample (with uncertainty between 1,2 and 4,9 cfu/25 g). It is possible to use a higher amount of Salmonella suspension for spiking: for example, if using 800 microliters, the reference value will be still valid (3,92 cfu/25 g sample) with better chance to obtain positive results, but the uncertainty of the spiking level will be slightly above the 5 cfu/sample (between 1,92 and 7,84 cfu/25 g).

To place an order or receive technical assistance

Order/Customer Service: SigmaAldrich.com/order Technical Service: SigmaAldrich.com/techservice

© 2023 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved. MilliporeSigma, the vibrant M, Millipore and Vitroids are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

MilliporeSigma is the U.S. and Canada Life Science business of Merck KGaA, Darmstadt, Germany,

MilliporeSigma 400 Summit Drive Burlington, MA 01803