

1.16730

Supelco®

sigmaaldrich.com

Reflectoquant®
Blank Strips for RQflex® 20
Leerstäbchen für RQflex® 20

50 Strips /
Stäbchen

Brief instructions for use:

1. Analyze the colored sample using a **test strip** from the Reflectoquant® test kit in question:
measurement value A
2. Repeat the analysis using a **blank strip**:
blank value B

Corrected result = A - B

In the event that this correction is not possible, decolorize the sample prior to the analysis using e.g. Polyvinylpolypyrrolidone Divergan® RS, Cat. No. 1.07302 (see also the Instruction for Use of the respective test kit).

Kurzgebrauchsanleitung:

1. Gefärbte Probe unter Verwendung eines **Test-stäbchens** aus dem betreffenden Reflectoquant® Testsatz analysieren:
Messwert A
2. Analyse unter Verwendung eines **Leerstäbchens** wiederholen:
Blindwert B

Korrigierter Messwert = A - B

Sollte diese Korrektur nicht möglich sein, Probe vor der Analyse entfärben, z. B. mit Polyvinylpolypyrrolidon Divergan® RS, Art. 1.07302 (s. auch Gebrauchsanweisung des betreffenden Testsatzes).

Please observe the RQflex® manual.



©2025

Store cool and dry • (15 - 25 °C) • Kühl und trocken lagern