

GenElute™-E Single Spin Checklist for Cleanup Kits

EC600 EC700 EC800

Prepare before starting

- Set the microcentrifuge to 1000 x g.

Column preparation

- Vortex Spin Column and place in a 2 mL tube for 5 minutes before continuing.



Loosen screw cap
of Spin Column.

OR



Punch a hole in the cap with the
GenElute™-E Single Spin Cap Puncher.

- Snap off bottom closure. Place Spin Column back into 2 mL tube.
- Centrifuge 1 minute at 1000 x g to collect column buffer.
- Place column in a 1.5 mL tube.

Purification of DNA

- Transfer 80 - 110 µL of sample to column.
- Centrifuge 1 minute at 1000 x g to collect nucleic acid.
- Collected nucleic acid is ready to use.

The life science business of Merck operates
as MilliporeSigma in the U.S. and Canada.

Merck, GenElute and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.
© 2022 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.

MERCK