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| ISO 10272による指定 | | FDA-BAM M30a による指定 | | GranuCult™ mCCDA基礎培地 (ISO 10272準拠) | |
|-----------------------------------|-----------|-----------------------------------|-----------|---------------------------------------|-----------|
| Meat Extract | 10 g/l | Meat Extract | 10 g/l | Meat Extract | 10 g/l |
| Enzymatic Digest of Animal Tissue | 10 g/l | Enzymatic Digest of Animal Tissue | 10 g/l | Enzymatic Digest of Animal Tissue | 10 g/l |
| Enzymatic Digest of Casein | 3 g/l | Enzymatic Digest of Casein | 3 g/l | Enzymatic Digest of Casein | 3 g/l |
| Charcoal | 4 g/l | Charcoal | 4 g/l | Charcoal | 4 g/l |
| Sodium Chloride | 5 g/l | Sodium Chloride | 5 g/l | Sodium Chloride | 5 g/l |
| Sodium Deoxycholate | 1 g/l | Sodium Deoxycholate | 1 g/l | Sodium Deoxycholate | 1 g/l |
| Sodium Pyruvate | 0.25 g/l | Sodium Pyruvate | 0.25 g/l | Sodium Pyruvate | 0.25 g/l |
| Iron (II) Sulfate | 0.25 g/l | Iron (II) Sulfate | 0.25 g/l | Iron (II) Sulfate | 0.25 g/l |
| - | - | Yeast Extract* | 2 g/l | - | - |
| Agar | 8-18 g/l | Agar | 8-18 g/l | Agar-Agar** | 12 g/l |
| 水 | 1000 ml/l | 水 | 1000 ml/l | 水 | n/a |
| 25°CのpH | 7.4 ± 0.2 | 25°CのpH | 7.4 ± 0.2 | 25°CのpH | 7.4 ± 0.2 |
| オートクレーブ滅菌後に添加されるサプリメント | | | | | |
| Cefoperazone | 0.032 g/l | | | Cefoperazone | 0.032 g/l |
| Amphotericin B | 0.01 g/l | | | Amphotericin B | 0.01 g/l |

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Merck Millipore, the M logo, GranuCult, Anaeroclip, Anaerocult and Singlepath are registered trademarks of Merck KGaA, Darmstadt, Germany. ATCC is a registered trademark of ATCC, Manassas, VA, USA. Lit. No. TN100070JA00

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Bolton, F.J. and Robertson, L. (1982): A selective medium for isolating *Campylobacter jejuni/coli*. J. Clin. Path. **35**: 462-467.

Bolton, F.J., Coates, D., Hinchcliffe, P.M. and Robertson, L. (1983): Comparison of selective media for isolation of *Campylobacter jejuni/coli*. J. Clin. Path. **36**: 78-83.

Bolton, F.J., Hutchinson, D.N. and Coates, D. (1984): Blood-free selective medium for isolation of *Campylobacter jejuni* from feces. J. Clin. Microbiol. **19**, 169-171.

Corry J.E.L. and Atabay, H.I. (2012): Culture media for the isolation of *Campylobacters*, *Helicobacters* and *Arcobacters*. In: Handbook of Culture Media for Food and Water Microbiology. (Corry, J.E.L., Curtis, G.D.W. and Baird, R.M. eds), pp. 403-435. Royal Society of Chemistry, Cambridge, UK.

FDA-BAM (2001): Chapter No. 7: *Campylobacter*. U.S. Food and Drug Administration – Bacteriological Analytical Manual.

ISO International Standardisation Organisation. Microbiology of food, animal feed and water – Horizontal method for detection and enumeration of *Campylobacter spp.* – Part 1: Detection method. EN ISO 10272-1:2006.

ISO International Standardisation Organisation. Microbiology of food and animal feeding stuffs – Horizontal method for detection and enumeration of *Campylobacter spp.* – Part 2: Colony-count technique EN ISO/TS 10272:2006.

ISO International Standardisation Organisation. Microbiology of food, animal feed and water – Preparation, production, storage and performance testing of culture media. EN ISO 11133:2014.

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