

## Product Information

# HRV-3C Protease

*N*-Terminal His-tagged recombinant protein

**SAE0045**

## Product Description

Synonym: Rhinovirus (serotype 14 LP) genomic region encoding protease

HRV-3C protease from human Rhinovirus Type 14 is a protease that specifically cleaves within the following eight-residue recognition sequence:

↓

Leu-Glu-Val-Leu-Phe-Gln-Gly-Pro

Proteolytic cleavage occurs between the Gln and Gly residues.<sup>1</sup> HRV-3C protease is a useful tool to cleave recombinant proteins that are expressed as fusion proteins with this sequence between the carrier domain and the protein of interest.<sup>2</sup>

This recombinant HRV-3C protease has a molecular mass of ~21 kDa. It contains a six-histidine tag (His-tag) to allow for removal by IMAC chromatography, such as with HIS-Select® or His•Bind® products. Several theses<sup>3-4</sup> and references<sup>5-6</sup> cite use of this SAE0045 HRV-3C protease product in their research protocols.<sup>3-4</sup>

## Reagent

The product is supplied in aqueous buffer, in a protein concentration range of 0.8-1.2 mg/mL, containing 50 mM Trizma-HCl (pH 7.5), 0.15 M NaCl, 1 mM TCEP, and 50% (v/v) glycerol. For the exact concentration, please refer to the lot-specific Certificate of Analysis for your specific lot number.

One unit of HRV-3C protease is defined as the amount of enzyme needed to digest 1 nmole of the substrate peptide H-Glu-Ala-Leu-Phe-Gln-pNA per hour at 0 °C, in a reaction buffer containing 25 mM HEPES (pH 7.5), 150 mM NaCl, 1 mM EDTA, and 1 mM DTT.

## Storage/Stability

The product retains activity for at least 2 years when stored at -20 °C.

## Procedure

HRV-3C protease is active under a wide range of pH values, ionic strengths, and temperatures. This protease retains high activity even at 0 °C, making it an optimal choice for temperature-sensitive proteins. However, the activity toward substrate proteins may differ depending on the substrate identity and reaction conditions.

The presence of low concentrations of a reducing agent in the reaction buffer is highly desirable, to keep the enzyme active in prolonged incubations. It is recommended to use 0.2-1 mM DTT in the reaction buffer for optimal results.

A good starting point for optimization is to use 1 µg of HRV-3C protease per 100 µg of target protein for 1 hour at 0-8 °C, or 1 µg of HRV-3C protease per 500 µg at 0-8 °C for 12-24 hours.

Temperatures up to 30 °C can be used for faster digestion. Protease activity is ~5× higher at 30 °C versus 0-8 °C. However, protease and substrate stability might be compromised.

## Precautions and Disclaimer

For R&D use only. Not for drug, household, or other uses. Please consult the Safety Data Sheet for information regarding hazards and safe handling practices.

## References

1. Cordingley, M.G. *et al.*, *J. Biol. Chem.*, **265**(16), 9062-9065 (1990).
2. Waugh, D.S., *Protein Expr. Purif.*, **80**(2), 283-293 (2011).
3. Aguilera, Marcos Moreno, "Estudio de Nrn1 en la patología de esquizofrenia" ("Study of Nrn1 in the pathology of schizophrenia"). Universidad de Barcelona, M.Sc. thesis, p. 9 (2017).
4. Gao, Mengjin, "The expression and purification of zinc metalloprotease STE24 (ZMPSTE24) and study of the interactions between the protein-inhibitor binding". McGill University, M.Sc. thesis, p. 41 (2021).
5. Lindesmith, L.C. *et al.*, *Immunity*, **50**(6), 1530-1541.e8 (2019).
6. Yukawa, M. *et al.*, *J. Exp. Med.*, **217**(1), e20182009 (2020).

## Notice

We provide information and advice to our customers on application technologies and regulatory matters to the best of our knowledge and ability, but without obligation or liability. Existing laws and regulations are to be observed in all cases by our customers. This also applies in respect to any rights of third parties. Our information and advice do not relieve our customers of their own responsibility for checking the suitability of our products for the envisaged purpose.

The information in this document is subject to change without notice and should not be construed as a commitment by the manufacturing or selling entity, or an affiliate. We assume no responsibility for any errors that may appear in this document.

### Technical Assistance

Visit the tech service page at [SigmaAldrich.com/techservice](https://SigmaAldrich.com/techservice).

### Standard Warranty

The applicable warranty for the products listed in this publication may be found at [SigmaAldrich.com/terms](https://SigmaAldrich.com/terms).

### Contact Information

For the location of the office nearest you, go to [SigmaAldrich.com/offices](https://SigmaAldrich.com/offices).

The life science business of Merck operates  
as MilliporeSigma in the U.S. and Canada.

Merck and Sigma-Aldrich are trademarks of Merck KGaA, Darmstadt, Germany or its affiliates. All other trademarks are the property of their respective owners. Detailed information on trademarks is available via publicly accessible resources.

© 2021 Merck KGaA, Darmstadt, Germany and/or its affiliates. All Rights Reserved.  
SAE0045pis Rev 01/22 DT,GCY,SM,MAM

