

## Product Information

### Streptolysin O from *Streptococcus pyogenes*

Catalog Numbers **S5265** and **S0149**

Storage Temperature 2–8 °C

CAS RN 98072-47-0

Synonym: SLO

#### Product Description

Streptolysin O (SLO) is a single polypeptide chain that is not glycosylated. It has a molecular mass of ~69 kDa<sup>1,2</sup> and an isoelectric point (pI) of 6.0-6.4.<sup>1</sup>

Streptolysin O is an immunogenic, oxygen-labile toxin, which is reversibly activated by dithiothreitol.<sup>2</sup> It is released into the extracellular medium along with other toxins, including streptolysin S, during the growth of most strains of group A and many strains of groups C and G *Streptococci*.<sup>2,3</sup> Streptolysin S differs from SLO in that it is oxygen-stable, nonimmunogenic and only active when associated with a carrier protein.<sup>4</sup> SLO may be used for cell permeabilization or hemolysis. Erythrocytes from different animal species have significantly different susceptibility to hemolysis by SLO.<sup>2</sup> It has been used to introduce oligonucleotides into cultured eukaryotic cells by permeabilization.<sup>4</sup> SLO has also been used to permeabilize human T lymphocytes.<sup>5</sup>

Kinetic studies have been performed on partially purified SLO. Cholesterol was found to inhibit hemolytic activity when added prior to incubation of SLO with rabbit erythrocytes; however, no inhibition was observed when cholesterol was added after the initiation of hemolysis.<sup>6</sup>

The products are lyophilized from a 10 mM Tris-Cl buffer solution containing 3 mM sodium azide, 5 mM sodium EDTA, and 1 mM PMSF. The protein content is 3-9% and each vial contains less than 1 mg solid (25,000–50,000 units/vial). S0149 has been  $\gamma$ -irradiated.

Unit Definition: One unit will cause 50% lysis of 50  $\mu$ l of a 2% human red blood cell suspension in phosphate buffered saline, pH 7.4, at 37 °C for 30 minutes.

#### Precautions and Disclaimer

This product is for R&D use only, not for drug, household, or other uses. Please consult the Material Safety Data Sheet for information regarding hazards and safe handling practices.

#### Preparation Instructions

SLO is soluble in cold water (1 mg/ml). Reconstitute by adding water into the vial and gently rotating.

After reconstitution, the solution should be stored at –20 °C in aliquots and any unused portions should be discarded. SLO is readily oxidized in solution. The activity is significantly increased in the presence of reducing agents, 20 mM cysteine<sup>3</sup> or 10 mM dithiothreitol.<sup>1</sup> A SLO solution will lose ~50% activity within 10 days when stored at 2-8 °C. It is recommended to freeze solutions at a minimal concentration of 0.2 mg/ml.

#### Storage/Stability

It is recommended to store Streptolysin O at 2–8 °C. The lyophilized products retain activity for up to three years when stored properly.

#### References

1. Bhakdi, S. *et al.*, *Infection and Immunity*, **46**, 394 (1984).
2. Alouf, J.E., *Pharmacol. Ther.*, **11**, 661-717 (1980).
3. Alouf, J.E., and Geoffroy, C., *Methods in Enzymology*, **165**, 52 (1988).
4. Barry, E. *et al.*, *Biotechniques*, **15**, 1016 (1993).
5. Graves, J.D. *et al.*, *Biochem. J.*, **265**, 407-413 (1990).
6. Kanbayashi, Y. *et al.*, *J. Biochem.*, **71**, 227-237 (1972).

EM,VC,NDH,TA,MAM 04/08-1

Sigma brand products are sold through Sigma-Aldrich, Inc.

Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications. Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply. Please see reverse side of the invoice or packing slip.