

OGYE Selective Supplement

Additive for the preparation of OGYE Selective Agar.

Mode of Action

OGYE Selective Supplement contains oxytetracycline in lyophilized form. It suppresses the growth of accompanying bacterial flora during culturing yeasts and moulds.

Composition (per vial)

Oxytetracycline in a buffered base 0.05 g.

Experimental Procedure

The lyophilisate is dissolved in the original vial by adding 10 ml of sterile, distilled water.

In the preparation of OGYE Agar, the dissolved content of one vial is evenly mixed into 500 ml sterile, still liquid medium cooled to 45-50 °C.

Ordering Information

Product	Ordering No.	Pack size
OGYE Selective Supplement	1.09877.0010	1 x 10 vials